

Firmware Revision for 1.9.5.202306010915

Firmware Change Log:

- Implement Current Limit as a variable through firmware and API
- Mirror naming convention of Filament 1 and Filament 2 to Filament A and Filament B in firmware and API
- Add a button to degas ion source to KJLC WebUI
- Change units during leak check
- Fix: Graph rescales incorrectly

To update the firmware:

- Click the Gear Icon in the top right of the KJLC RGA Web UI
- Click the Green Firmware upload button
- Upload the LINXON.bin file first
- Upload the httproot.mar file second
- Clear the cache in your browser and power cycle your RGA